

# Арматура Cleanfit CPA875

## Техническая информация

По вопросам продаж и поддержки обращайтесь:

Алматы (7273)495-231  
Ангарск (3955)60-70-56  
Архангельск (8182)63-90-72  
Астрахань (8512)99-46-04  
Барнаул (3852)73-04-60  
Белгород (4722)40-23-64  
Благовещенск (4162)22-76-07  
Брянск (4832)59-03-52  
Владивосток (423)249-28-31  
Владикавказ (8672)28-90-48  
Владимир (4922)49-43-18  
Волгоград (844)278-03-48  
Вологда (8172)26-41-59  
Воронеж (473)204-51-73  
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06  
Ижевск (3412)26-03-58  
Иркутск (395)279-98-46  
Казань (843)206-01-48  
Калининград (4012)72-03-81  
Калуга (4842)92-23-67  
Кемерово (3842)65-04-62  
Киров (8332)68-02-04  
Коломна (4966)23-41-49  
Кострома (4942)77-07-48  
Краснодар (861)203-40-90  
Красноярск (391)204-63-61  
Курск (4712)77-13-04  
Курган (3522)50-90-47  
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13  
Москва (495)268-04-70  
Мурманск (8152)59-64-93  
Набережные Челны (8552)20-53-41  
Нижний Новгород (831)429-08-12  
Новокузнецк (3843)20-46-81  
Ноябрьск (3496)41-32-12  
Новосибирск (383)227-86-73  
Омск (3812)21-46-40  
Орел (4862)44-53-42  
Оренбург (3532)37-68-04  
Пенза (8412)22-31-16  
Петрозаводск (8142)55-98-37  
Псков (8112)59-10-37  
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15  
Рязань (4912)46-61-64  
Самара (846)206-03-16  
Санкт-Петербург (812)309-46-40  
Саратов (845)249-38-78  
Севастополь (8692)22-31-93  
Саранск (8342)22-96-24  
Симферополь (3652)67-13-56  
Смоленск (4812)29-41-54  
Сочи (862)225-72-31  
Ставрополь (8652)20-65-13  
Сургут (3462)77-98-35  
Сыктывкар (8212)25-95-17  
Тамбов (4752)50-40-97  
Тверь (4822)63-31-35

Тольятти (8482)63-91-07  
Томск (3822)98-41-53  
Тула (4872)33-79-87  
Тюмень (3452)66-21-18  
Ульяновск (8422)24-23-59  
Улан-Удэ (3012)59-97-51  
Уфа (347)229-48-12  
Хабаровск (4212)92-98-04  
Чебоксары (8352)28-53-07  
Челябинск (351)202-03-61  
Череповец (8202)49-02-64  
Чита (3022)38-34-83  
Якутск (4112)23-90-97  
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

эл.почта: [ehr@nt-rt.ru](mailto:ehr@nt-rt.ru) || сайт: <https://endcounters.nt-rt.ru/>

## Manual or automatic retractable assembly Cleanfit CPA875

Sterilizable assembly for the life sciences and food & beverage industries



### Benefits:

- Highest process safety: The patented design of the dynamic sealing prevents contamination.
- Simple operation: Remove, sterilize and calibrate the sensor while the process is running – aseptic and hygienically safe.
- No cross-contamination: Sterile design according to EHEDG and ASME BPE as well as sealings according to FDA and USP class VI.
- Flexible: Individual adaptation to your process requirements based on variety of process connections.
- Optional double service chamber: Sealing between sterile and non-sterile part moves dynamically with the sensor guide – just like in a syringe. The sterile space and the process medium remain untouched.

### Specs at a glance

- **Process temperature** -10 to 140 °C (14 to 284 °F)
- **Process pressure** Pneumatic drive: 16 bar up to 140 °C (232 psi up to 284 °F) Manual drive: 8 bar up to 140 °C (116 psi up to 284 °F)

**Field of application:** Completely sterilizable, the Cleanfit CPA875 retractable assembly guarantees maximum protection from contamination making it ideal for use in life science and food & beverage applications. Fully certified (EHEDG and ASME BPE) and made of FDA-recommended materials, the CPA875 is the right choice where hygiene counts!

### Features and specifications

## pH

**Measuring principle**

Potentiometric

---

**Application**

Food and beverages, biotechnology, life sciences, special chemicals

---

**Installation**

Retractable process assembly

---

**Characteristic**

Open and closed tanks, piping

---

**Design**

Sterile and hygienic stile

Manual or pneumatic

---

**Material**

Holder: stainless steel 1.4435

Seals: EPDM-FDA, FKM-FDA, FFKM-FDA

---

**Dimension**

Immersion depth: 16.1 to 81.4 mm (0.63 to 3.20 inch), depending on process adaption

---

**Process temperature**

-10 to 140 °C (14 to 284 °F)

---

**Process pressure**

Pneumatic drive: 16 bar up to 140 °C (232 psi up to 284 °F)

Manual drive: 8 bar up to 140 °C (116 psi up to 284 °F)

---

**Connection**

Aseptic DN25 clampable, Aseptic DN50 screw-in, Clamp ISO 2852, ASME BPE, Varivent N

---

## Oxygen

**Measuring principle**

Sensor ORP / Redox

---

## Oxygen

**Application**

Food and beverages, biotechnology, life sciences, special chemicals

---

**Installation**

Retractable process assembly

---

**Characteristic**

Open and closed tanks, piping

---

**Design**

Sterile and hygienic stile

Manual or pneumatic

---

**Material**

Holder: stainless steel 1.4435

Seals: EPDM-FDA, FKM-FDA, FFKM-FDA

---

**Dimension**

Immersion depth: 16.1 to 81.4 mm (0.63 to 3.20 inch), depending on process adaption

---

**Process temperature**

-10 to 140 °C (14 to 284 °F)

---

**Process pressure**

Pneumatic drive: 16 bar up to 140 °C (232 psi up to 284 °F)

Manual drive: 8 bar up to 140 °C (116 psi up to 284 °F)

---

**Connection**

Aseptic DN25 clampable, Aseptic DN50 screw-in, Clamp ISO 2852, ASME BPE, Varivent N

---

## По вопросам продаж и поддержки обращайтесь:

Алматы (7273)495-231  
Ангарск (3955)60-70-56  
Архангельск (8182)63-90-72  
Астрахань (8512)99-46-04  
Барнаул (3852)73-04-60  
Белгород (4722)40-23-64  
Благовещенск (4162)22-76-07  
Брянск (4832)59-03-52  
Владивосток (423)249-28-31  
Владикавказ (8672)28-90-48  
Владимир (4922)49-43-18  
Волгоград (844)278-03-48  
Вологда (8172)26-41-59  
Воронеж (473)204-51-73  
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06  
Ижевск (3412)26-03-58  
Иркутск (395)279-98-46  
Казань (843)206-01-48  
Калининград (4012)72-03-81  
Калуга (4842)92-23-67  
Кемерово (3842)65-04-62  
Киров (8332)68-02-04  
Коломна (4966)23-41-49  
Кострома (4942)77-07-48  
Краснодар (861)203-40-90  
Красноярск (391)204-63-61  
Курск (4712)77-13-04  
Курган (3522)50-90-47  
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13  
Москва (495)268-04-70  
Мурманск (8152)59-64-93  
Набережные Челны (8552)20-53-41  
Нижний Новгород (831)429-08-12  
Новокузнецк (3843)20-46-81  
Ноябрьск (3496)41-32-12  
Новосибирск (383)227-86-73  
Омск (3812)21-46-40  
Орел (4862)44-53-42  
Оренбург (3532)37-68-04  
Пенза (8412)22-31-16  
Петрозаводск (8142)55-98-37  
Псков (8112)59-10-37  
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15  
Рязань (4912)46-61-64  
Самара (846)206-03-16  
Санкт-Петербург (812)309-46-40  
Саратов (845)249-38-78  
Севастополь (8692)22-31-93  
Саранск (8342)22-96-24  
Симферополь (3652)67-13-56  
Смоленск (4812)29-41-54  
Сочи (862)225-72-31  
Ставрополь (8652)20-65-13  
Сургут (3462)77-98-35  
Сыктывкар (8212)25-95-17  
Тамбов (4752)50-40-97  
Тверь (4822)63-31-35

Тольятти (8482)63-91-07  
Томск (3822)98-41-53  
Тула (4872)33-79-87  
Тюмень (3452)66-21-18  
Ульяновск (8422)24-23-59  
Улан-Удэ (3012)59-97-51  
Уфа (347)229-48-12  
Хабаровск (4212)92-98-04  
Чебоксары (8352)28-53-07  
Челябинск (351)202-03-61  
Череповец (8202)49-02-64  
Чита (3022)38-34-83  
Якутск (4112)23-90-97  
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

эл.почта: [ehr@nt-rt.ru](mailto:ehr@nt-rt.ru) || сайт: <https://endcounters.nt-rt.ru/>