

# Датчики рН, ОВП Memosens CPS171D

## Техническая информация

По вопросам продаж и поддержки обращайтесь:

Алматы (7273)495-231  
Ангарск (3955)60-70-56  
Архангельск (8182)63-90-72  
Астрахань (8512)99-46-04  
Барнаул (3852)73-04-60  
Белгород (4722)40-23-64  
Брянск (4832)59-03-52  
Владивосток (423)249-28-31  
Владикавказ (8672)28-90-48  
Владимир (4922)49-43-18  
Волгоград (844)278-03-48  
Вологда (8172)26-41-59  
Воронеж (473)204-51-73  
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06  
Ижевск (3412)26-03-58  
Иркутск (395)279-98-46  
Казань (843)206-01-48  
Калининград (4012)72-03-81  
Калуга (4842)92-23-67  
Кемерово (3842)65-04-62  
Киров (8332)68-02-04  
Коломна (4966)23-41-49  
Кострома (4942)77-07-48  
Краснодар (861)203-40-90  
Красноярск (391)204-63-61  
Курск (4712)77-13-04  
Курган (3522)50-90-47  
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13  
Москва (495)268-04-70  
Мурманск (8152)59-64-93  
Набережные Челны (8552)20-53-41  
Нижний Новгород (831)429-08-12  
Новокузнецк (3843)20-46-81  
Ноябрьск (3496)41-32-12  
Новосибирск (383)227-86-73  
Омск (3812)21-46-40  
Орел (4862)44-53-42  
Оренбург (3532)37-68-04  
Пенза (8412)22-31-16  
Петрозаводск (8142)55-98-37  
Псков (8112)59-10-37  
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15  
Рязань (4912)46-61-64  
Самара (846)206-03-16  
Санкт-Петербург (812)309-46-40  
Саратов (845)249-38-78  
Севастополь (8692)22-31-93  
Саранск (8342)22-96-24  
Симферополь (3652)67-13-56  
Смоленск (4812)29-41-54  
Сочи (862)225-72-31  
Ставрополь (8652)20-65-13  
Сургут (3462)77-98-35  
Сыктывкар (8212)25-95-17  
Тамбов (4752)50-40-97  
Тверь (4822)63-31-35

Тольятти (8482)63-91-07  
Томск (3822)98-41-53  
Тула (4872)33-79-87  
Тюмень (3452)66-21-18  
Ульяновск (8422)24-23-59  
Улан-Удэ (3012)59-97-51  
Уфа (347)229-48-12  
Хабаровск (4212)92-98-04  
Чебоксары (8352)28-53-07  
Челябинск (351)202-03-61  
Череповец (8202)49-02-64  
Чита (3022)38-34-83  
Якутск (4112)23-90-97  
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

эл.почта: [ehr@nt-rt.ru](mailto:ehr@nt-rt.ru) || сайт: <https://endcounters.nt-rt.ru/>

# Digital pH sensor Memosens CPS171D

## Memosens glass electrode for bioreactors in the biotech industry and for the food industry



### Benefits:

- The sensor's exceptional accuracy, reproducibility and reliability helps you to keep the pH value in the optimum range for a maximized product yield.
- A specialized glass membrane and reference system makes the sensor CIP/SIP and autoclaving resistant (up to 140°C/284°F) offering maximum long-term stability.
- Thanks to the Memosens technology, the cable connection cannot be affected by moisture or corrosion: you can count on the sensor's highest availability and your batch is safe.
- The sensor protects your product quality with its certified biocompatibility according to USP class VI, FDA compliance of wetted parts, no cytotoxicity, bioreactivity or animal-based materials.
- The sensor provides you with process and sensor data that allows you to decide if the sensor is ready for the next batch or should be replaced - a big improvement for your product safety.
- Streamline maintenance of CPS171D with the Memobase Plus software that automatically stores and documents all relevant sensor and process data.
- Approved for hazardous area use.

### Specs at a glance

- **Measurement range** pH 0 to 14
- **Process temperature** 0 to 135 °C (32 to 275 °F)
- **Process pressure** 1 to 7 bar abs (14.5 to 101.5 psi)

**Field of application:** Keep the pH value of your batch in the optimum range and maximize your product yield with Memosens CPS171D. Thanks to its rugged design and long-term stability, it provides extremely

accurate, reproducible and reliable measuring values even after CIP/SIP and autoclaving. Memosens CPS171D is not the latest Memosens generation. To get information on the new Memosens CPS61E sensor with extended functionality, click [here](#).

## Features and specifications

pH

### Measuring principle

Potentiometric

### Application

Hygienic and sterile applications, bioreactor, fermenter, biotechnology, pharmaceutical industry, foods

### Characteristic

Digital pH electrodes for bioreactors and biotechnology production processes with ion trap for long-term stable reference

### Measurement range

pH 0 to 14

### Measuring principle

Ceramic diaphragm, ion-trap

### Design

All shaft lengths with temperature sensor  
Advanced gel technology  
Digital electrode with Memosens technology

### Material

Glass and ceramic

### Dimension

Diameter: 12 mm (0.47 inch)  
Shaft lengths: 120, 225, 360 and 425 mm  
(4.72, 8.86, 14.2 and 16.7 inch)

pH

**Process temperature**

0 to 135 °C (32 to 275 °F)

---

**Process pressure**

1 to 7 bar abs (14.5 to 101.5 psi)

---

**Temperature sensor**

NTC 30k

---

**Ex certification**

IECEX, ATEX, NEPSI, FM, CSA

---

**Connection**

Inductive, digital connection head with Memosens technology

---

## По вопросам продаж и поддержки обращайтесь:

Алматы (7273)495-231  
Ангарск (3955)60-70-56  
Архангельск (8182)63-90-72  
Астрахань (8512)99-46-04  
Барнаул (3852)73-04-60  
Белгород (4722)40-23-64  
Брянск (4832)59-03-52  
Владивосток (423)249-28-31  
Владикавказ (8672)28-90-48  
Владимир (4922)49-43-18  
Волгоград (844)278-03-48  
Вологда (8172)26-41-59  
Воронеж (473)204-51-73  
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06  
Ижевск (3412)26-03-58  
Иркутск (395)279-98-46  
Казань (843)206-01-48  
Калининград (4012)72-03-81  
Калуга (4842)92-23-67  
Кемерово (3842)65-04-62  
Киров (8332)68-02-04  
Коломна (4966)23-41-49  
Кострома (4942)77-07-48  
Краснодар (861)203-40-90  
Красноярск (391)204-63-61  
Курск (4712)77-13-04  
Курган (3522)50-90-47  
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13  
Москва (495)268-04-70  
Мурманск (8152)59-64-93  
Набережные Челны (8552)20-53-41  
Нижний Новгород (831)429-08-12  
Новокузнецк (3843)20-46-81  
Ноябрьск (3496)41-32-12  
Новосибирск (383)227-86-73  
Омск (3812)21-46-40  
Орел (4862)44-53-42  
Оренбург (3532)37-68-04  
Пенза (8412)22-31-16  
Петрозаводск (8142)55-98-37  
Псков (8112)59-10-37  
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15  
Рязань (4912)46-61-64  
Самара (846)206-03-16  
Санкт-Петербург (812)309-46-40  
Саратов (845)249-38-78  
Севастополь (8692)22-31-93  
Саранск (8342)22-96-24  
Симферополь (3652)67-13-56  
Смоленск (4812)29-41-54  
Сочи (862)225-72-31  
Ставрополь (8652)20-65-13  
Сургут (3462)77-98-35  
Сыктывкар (8212)25-95-17  
Тамбов (4752)50-40-97  
Тверь (4822)63-31-35

Тольятти (8482)63-91-07  
Томск (3822)98-41-53  
Тула (4872)33-79-87  
Тюмень (3452)66-21-18  
Ульяновск (8422)24-23-59  
Улан-Удэ (3012)59-97-51  
Уфа (347)229-48-12  
Хабаровск (4212)92-98-04  
Чебоксары (8352)28-53-07  
Челябинск (351)202-03-61  
Череповец (8202)49-02-64  
Чита (3022)38-34-83  
Якутск (4112)23-90-97  
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(7172)727-132

Киргизия +996(312)96-26-47

эл.почта: [ehr@nt-rt.ru](mailto:ehr@nt-rt.ru) || сайт: <https://endcounters.nt-rt.ru/>